

**USING PROFILOMETRY  
FOR FAST 2D MEASUREMENT**



Prepared by  
**Andrea Novitsky**

## INTRODUCTION:


Whether the end application is production, quality control, or research and development, access to quick and precise 2D measurement is needed. Although 3D area measurement may ultimately also be used having both provides time saving and broadens instrument use. The Nanovea Non-Contact Profilometer is capable of 3D & 2D from the nano to macro scale providing broad range usability.

### IMPORTANCE OF PROFILOMETRY FOR FAST 2D MEASUREMENT

Unlike other techniques such as touch probes or interferometry, the Non Contact Profilometer, using axial chromatism, can measure nearly any surface, sample sizes can vary widely due to open staging and there is no sample preparation needed. Nano through macro range is obtained during surface measurement with zero influence from sample reflectivity or absorption, has advanced ability to measure high surface angles and there is no software manipulation of results. Easily measure any material: transparent, opaque, specular, diffusive, polished, rough etc. The technique of the Non Contact Profilometer provides an ideal, broad and user friendly capability to maximize surface studies when with the benefits of combined 2D & 3D capability.


### MEASUREMENT OBJECTIVE

In this application the Nanovea ST400 Profilometer is used to measure roughness, step heights and other dimensions on a machined part. All of these calculations are made from a single 2D measurement that takes less than 5 seconds to obtain. This speed and ease of use combined with highly accurate data makes the Nanovea Profilometer a valuable tool for a wide range of applications.

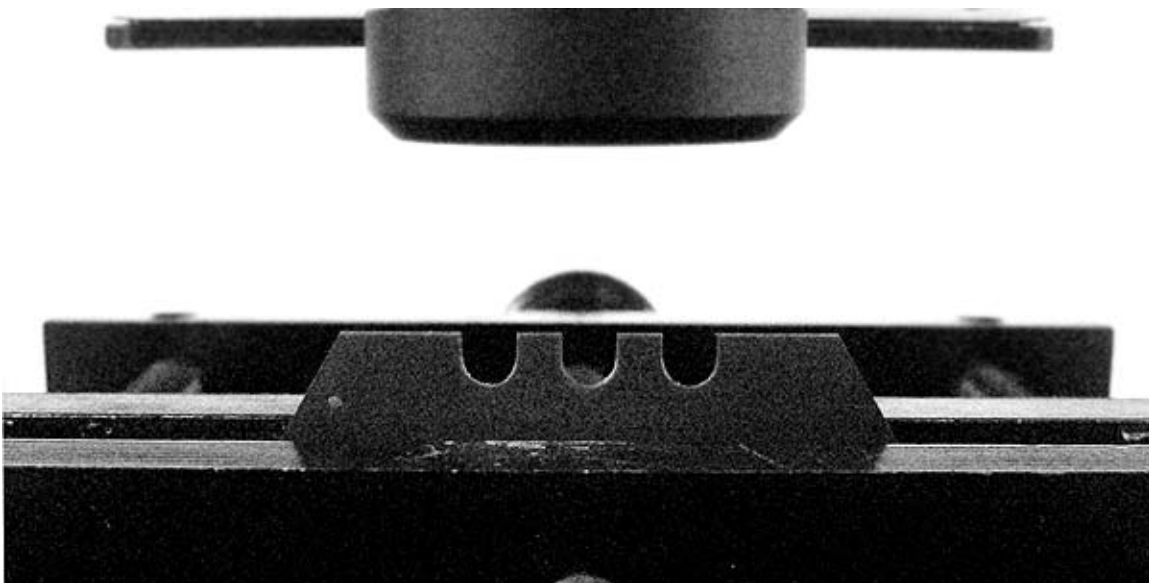

## RESULTS: 2D Measurement

The roughness, step height and contour were used to calculate various dimensions from a single line profile measurement.

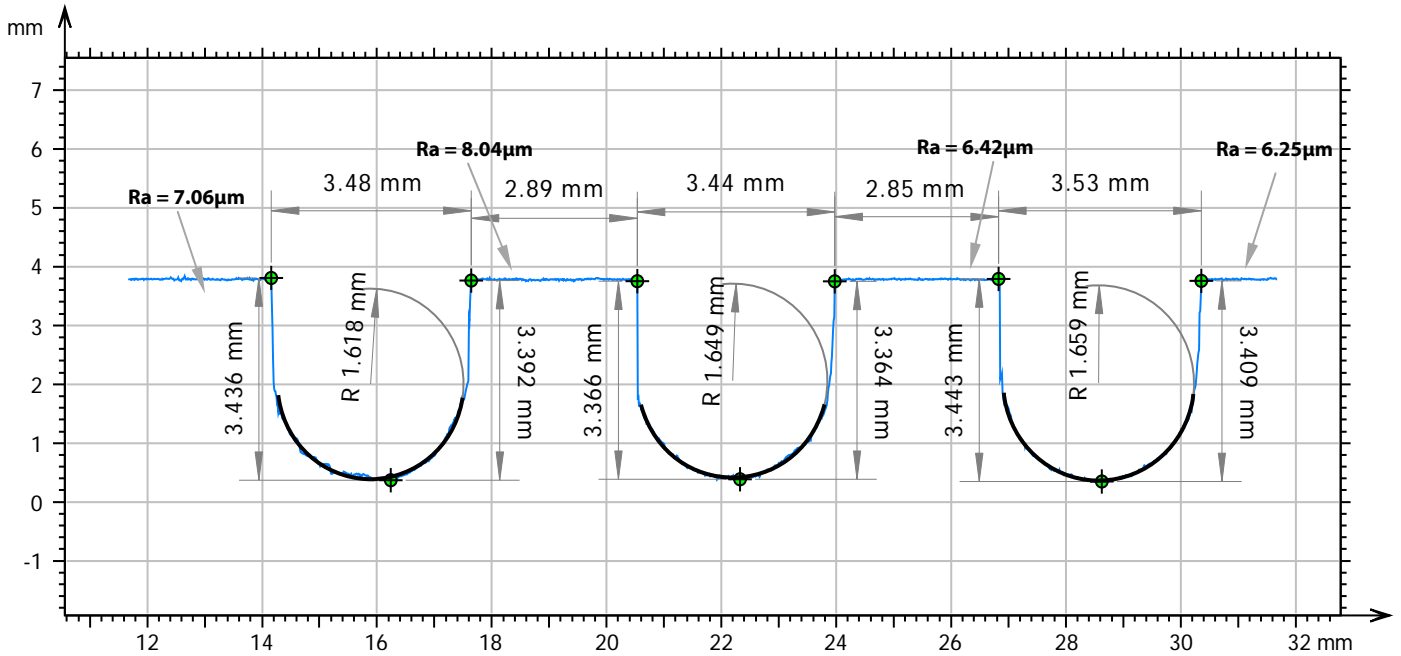


## CONCLUSION:

In this application, we have shown how the Nanovea Non-Contact Profilometer can precisely measure roughness, various heights, widths, and other dimensions from a single 2D measurement that takes seconds to obtain. The data shows an average step height of 3.418mm, an average radius of 1.642mm, as well as other dimensions. These measurements are relevant to a huge range of applications, from precision machined parts to biomedical products to electronics. The data shown here represents only a portion of the calculations available in the analysis software.

Learn more about the [Nanovea Profilometer](#) or [Lab Services](#)

## MEASUREMENT PRINCIPLE:

The Chromatic Confocal technique uses a white light source, where light passes through an objective lens with a high degree of chromatic aberration. The refractive index of the objective lens will vary in relation to the wavelength of the light. In effect, each separate wavelength of the incident white light will re-focus at a different distance from the lens (different height). When the measured sample is within the range of possible heights, a single monochromatic point will be focalized to form the image. Due to the confocal configuration of the system, only the focused wavelength will pass through the spatial filter with high efficiency, thus causing all other wavelengths to be out of focus. The spectral analysis is done using a diffraction grating. This technique deviates each wavelength at a different position, intercepting a line of CCD, which in turn indicates the position of the maximum intensity and allows direct correspondence to the Z height position.

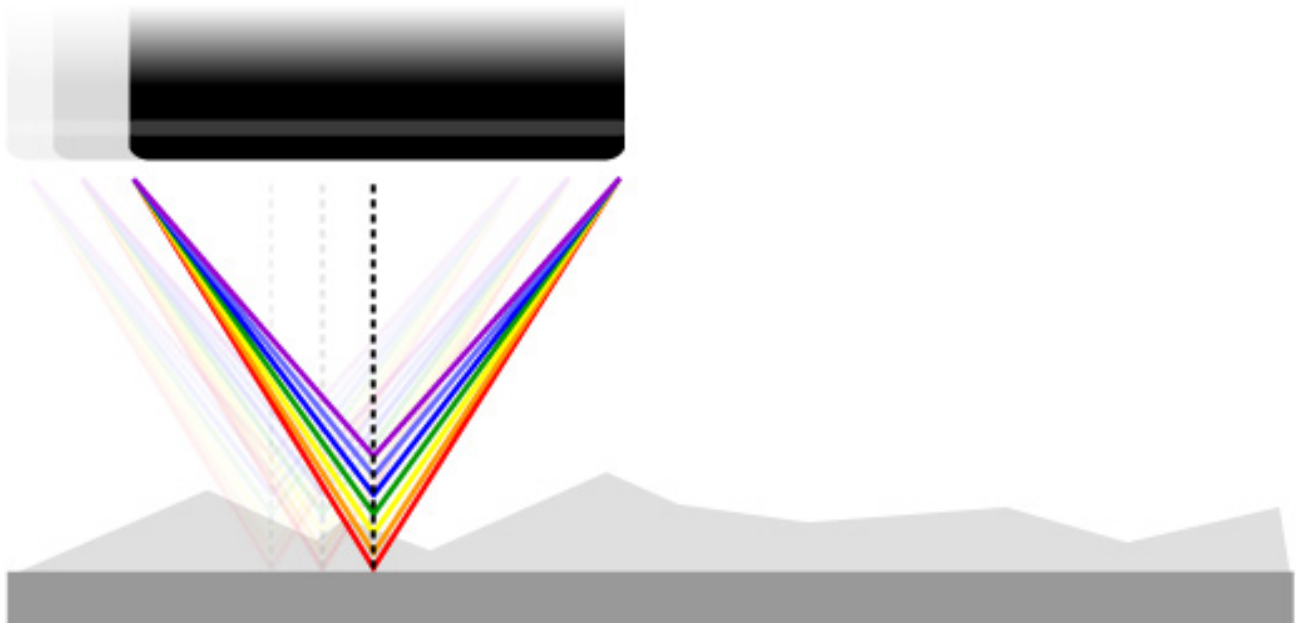

Unlike the errors caused by probe contact or the manipulative Interferometry technique, Chromatic Confocal technology measures height directly from the detection of the wavelength that hits the surface of the sample in focus. It is a direct measurement with no mathematical software manipulation. This provides unmatched accuracy on the surface measured because a data point is either measured accurately without software interpretation or not at all. The software completes the unmeasured point but the user is fully aware of it and can have confidence that there are no hidden artifacts created by software guessing.

Nanovea optical pens have zero influence from sample reflectivity or absorption. Variations require no sample preparation and have advanced ability to measure high surface angles. Capable of large Z measurement ranges. Measure any material: transparent or opaque, specular or diffusive, polished or rough. Measurement includes: Profile Dimension, Roughness Finish Texture, Shape Form Topography, Flatness Warpage Planarity, Volume Area, Step-Height Depth Thickness and many others.

## DEFINITION OF HEIGHT PARAMETERS

Height Parameter		Definition
Sa	Arithmetical Mean Height	Mean surface roughness. $Sa = \frac{1}{A} \iint_A  z(x, y)  dx dy$
Sq	Root Mean Square Height	Standard deviation of the height distribution, or RMS surface roughness. $Sq = \sqrt{\frac{1}{A} \iint_A z^2(x, y) dx dy}$ <p>Computes the standard deviation for the amplitudes of the surface (RMS).</p>
Sp	Maximum Peak Height	Height between the highest peak and the mean plane.
Sv	Maximum Pit Height	Depth between the mean plane and the deepest valley.
Sz	Maximum Height	Height between the highest peak and the deepest valley.
Ssk	Skewness	Skewness of the height distribution. $Ssk = \frac{1}{Sq^3} \left[ \frac{1}{A} \iint_A z^3(x, y) dx dy \right]$ <p>Skewness qualifies the symmetry of the height distribution. A negative Ssk indicates that the surface is composed of mainly one plateau and deep and fine valleys. In this case, the distribution is sloping to the top. A positive Ssk indicates a surface with a lot of peaks on a plane. Therefore, the distribution is sloping to the bottom.</p> <p>Due to the large exponent used, this parameter is very sensitive to the sampling and noise of the measurement.</p>
Sku	Kurtosis	Kurtosis of the height distribution. $Sku = \frac{1}{Sq^4} \left[ \frac{1}{A} \iint_A z^4(x, y) dx dy \right]$ <p>Kurtosis qualifies the flatness of the height distribution.</p> <p>Due to the large exponent used, this parameter is very sensitive to the sampling and noise of the measurement.</p>
Spar	Projected Area	Projected surface area.
Sdar	Developed Area	Developed surface area.